

University of Wisconsin–Madison
Chemistry Department

Varian NMR User's Guide

by: Charles G. Fry

(updated February 24, 2003)

Note: This guide provides an introduction to use of Varian equipment at the UWChemMRF. *This guide is not intended in any way to be a replacement to the excellent Varian documentation!* All students should refer regularly to the Varian VNMR Liquids Users Guide for learning and the Varian VNMR Command Reference Guide for specific guidance. All the Varian documentation is available in both hardcopy and on-line.

*Copyright © 1996-2003 Charles G. Fry
All Rights Reserved.*

Table of Contents


UW Chemistry Magnetic Resonance Facility (UWChemMRF)	4
I. Facility Layout (2 nd floor Matthews)	4
SGIs – NATOTH (Avance host computer) and GQUAN (for off-line data workup)	5
II. Facility Personnel	5
Quick Guide for New Users	6
I. Login	6
II. Setup	6
A. 1 st Time:	6
B. Parameters:	6
C. Variable Temperature	7
III. Locking and Shimming	7
A. Sample Prep:	7
B. Locking:	7
C. Shimming:	8
IV. Probe Tuning and ¹ H Calibration	8
V. Acquisition	8
VI. Saving and Deleting Data, and Backups	9
VII. Logging Out	9
VIII. Don'ts	9
Introduction to Unix and VNMR	10
I. Common Unix Commands	10
II. Common VNMR Commands, Parameters and Flags	12
III. VNMR Directory Structure	16
IV. Common Pulse Sequence Setup in VNMR	17
V. VNMR to Bruker-AM/AC Parameter Conversion Table	19
VI. Trouble-Shooting	25
Operation of Unity-500	27
I. Proper Exiting and Logging In and Initial VNMR Setup	27
II. Commands for First Time and Novice Users	27
III. Probe Changes	27
Table 1. Description of Probes on Unity-500 and Inova-500	29
Table 2. Calibrations of Probes on Unity-500	31
III. Probe Tuning	31
IV. Lock and shim	33
Table 3. Field and Lock Power Settings for Unity-500	33
Table 4. Major Shim Interactions on Unity-500	34
Table 5. Shim Sensitivities on Unity-500	34
Experiments on Unity-500	35
I. Normal 1d ¹ H Acquisition	35
D. ¹ H pw90 Calibration	36
E. 2 nd Order Shimming on a 500 MHz Spectrometer	37

F. 1d Data Workup and Plotting	37
II. X Acquisition (e.g., ^{31}P or ^{13}C)	39
Table 6. Common Decoupler Parameter Settings	40
D. ^{13}C (X) pw90 and $\{^1\text{H}\}$ Decoupler Calibrations	41
III. DEPT – Distortionless Enhancement by Polarization Transfer	43
IV. INEPT – Insensitive Nuclei Enhanced by Polarization Transfer	46
V. 1d NOE-Difference Spectroscopy	48
VI. Homonuclear 1d Decoupling	50
VII. COSY – 2d Homonuclear Correlation	51
E. 2d Data Workup and Plotting	52
VIII. DQ-COSY – Double-Quantum Filtered COSY (Phase-Sensitive)	53
IX. TOCSY (or HOHAHA) – Total Correlation Spectroscopy	55
X. Other COSY Variants	57
XI. NOESY – 2d NOE and Exchange Spectroscopy	58
XII. ROESY – 2d NOE Spec.; Mixing via a Spinlock in the Rotating Frame	60
XIII. HETCOR – 2d Heterocorrelation Spec. (“normal” X nucleus detection)	62
XIV. HMQC and HSQC – 2d Heteronuclear Spectroscopy (“inverse” experiment with ^1H detection)	63
XV. HMBC – 2d Long-Range Heterocorrelation Spec. (inverse detection)	65
XVI. Other Heteronuclear Correlation Experiments	66
XVII. Fitting Partially Overlapping Peaks – Deconvolution in VNMR	67
XIX. ^1H Spin-Lattice Relaxation, T_1	68
II. Rapid Determination of T_1 by Inversion-Recovery Null Method	68
III. Quantitative measurement of T_1 by Inversion Recovery Method	69
A. Comments	69
B. Acquisition Set-up	69
C. T_1 Analysis (see Varian Subject manual, Adv. 1d section)	69
XIX. Cross-Polarization/Magic-Angle-Spinning NMR	72
A. Preparation of MAS spectra	73
B. Simulation of spinning sideband patterns	74
C. Least-squares fitting of spinning sideband patterns	74
Data Backups from UNIX Workstations	76
I. Backups to Colorado Systems Tape Drive on Windows NT Network	76
A. Prepare tar file on UNIX workstations	76
B. Move file to NT server	76
C. Backup file to Colorado Systems tape drive	76
D. Recovering backup'ed up data	76
II. Backups to IOmega Zip Drive on Windows NT Network (currently installed on Morder, Terminus, Drazi)	76
A. Prepare a tar file	76
B. Backup file to IOmega Zip drive	77
D. Recovering backup'ed up data	77
III. Backups to Pinnacle Optical Disk (on Tango on 5 th floor)	77
A. Format a New Optical Disk	77
B. Create File System on New Optical Disk	77
C. Backup to Optical Disk	77
IV. Backups to Pinnacle CD Writer (on Twiddle on 9 th floor)	78

A. Transfer Data to Twiddle	78
B. Transfer Directories/Files to Easy-CD Window	78
Administration of Varian Spectrometers.....	80
I. Rebooting Sun's on Spectrometers (requires root privileges).....	80
A. Shutting down UNIX:	80
B. Powering Up UNIX:	80
Pulsed-Field Gradient Shimming With VNMR.....	81
I. General Discussion	81
II. Normal PFG Shimming Procedure.....	82
III. First-Time Use and Detailed Explanation of Parameters	83
A. Critical Parameters	84
B. Details of Correct Parameter Setup	85

UW Chemistry Magnetic Resonance Facility (UWChemMRF)

I. Facility Layout (2nd floor Matthews)

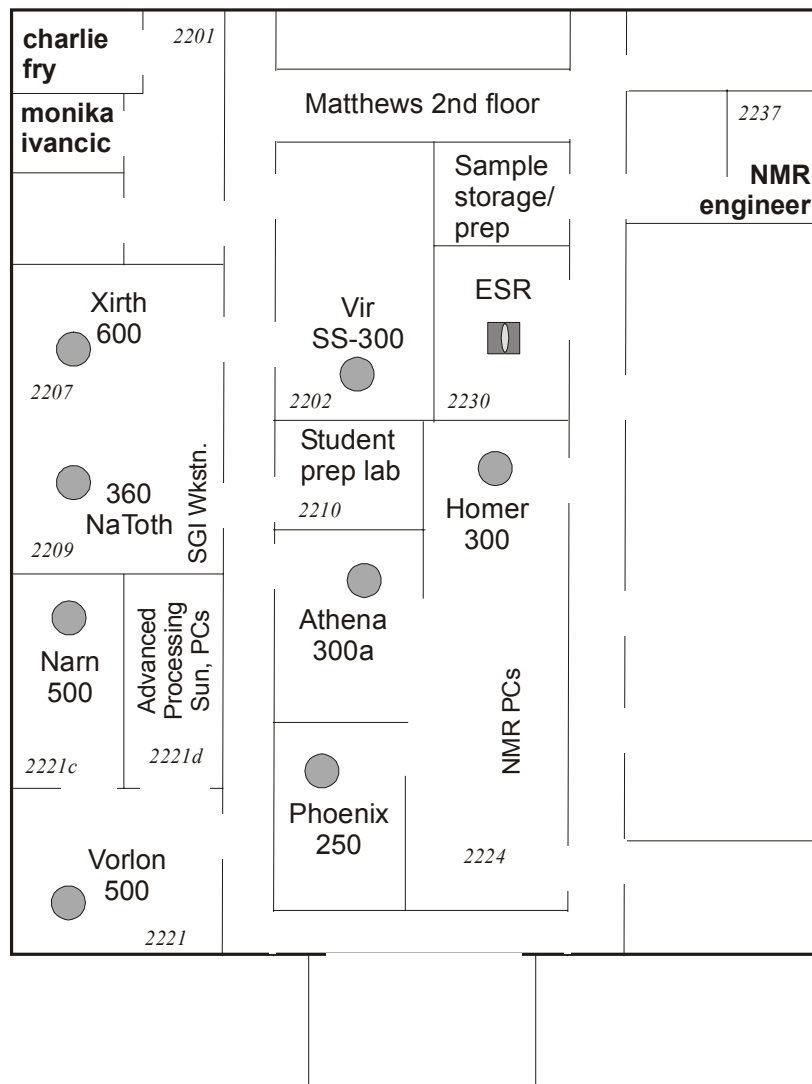

As of Feb. 2003

ATHENA	–	AC+ 300	routine $^1\text{H}/^{19}\text{F}/^{31}\text{P}/^{13}\text{C}$	–	sample changer, quad-nucleus probe
HOMER	–	AC+ 300	routine $^1\text{H}/^{13}\text{C}$	–	$^1\text{H}/^{13}\text{C}$ dedicated
PHOENIX	–	AC+ 250	routine BB VT	–	routine BB ($^{29}\text{Si}/^{11}\text{B}/^2\text{H}/^{199}\text{Hg}/\text{etc.}$), variable temperature
VIR	–	UNITY-300	solid-state NMR	–	conformational, motions, solid-state packing, catalysts, amorphous and glassy compounds
NATOTH	–	Avance-360	non-routine BB VT	–	long-term VT, kinetics, concentration limited samples; 5 and 10 mm BB probes, 5 mm inverse probe

NARN	–	UNITY-500 non-routine ^1H /BB VT	–	high-sensitivity, sample-limited ($^1\text{H} < 5 \text{ mg}$, $^{13}\text{C} < 15 \text{ mg}$), short-run, sophisticated experiments (e.g., HMQC, DQCOSY, gCOSY, gNOESY); limited access
VORLON	–	INOVA-500 inverse exps, 2D studies	–	long-term, sophisticated, gradient-enhanced experiments; combi-chem MAS probe; limited access
XIRTH	–	INOVA-600 long-term 2D studies	–	long-term, most sophisticated, gradient-enhanced experiments (e.g., NOESY/ROESY, HMQC, DQCOSY); limited access
ESR	–	ESP-300 electron spin resonance	–	paramagnetism, free-radical chemistry
PC's	–	Six PC's are located in rm. 2224 for use in working up data. The PCs connect to a Win2000 server, which uploads data once per minute to a SNAP server, nmrsnap.chem.wisc.edu (128.104.70.79). The SNAP server can be mapped to PCs or Macs within the department, or accessed via IE or Netscape. See the facility director for more information		
	–	All PCs have Acorn Software's NUTS 1D, H. Reich's WinDNMR, FTP and Zip software loaded. Some PCs have NUTS 2D Pro, ACDs HNMR and CNMR, gNMR, and Mestre-C software loaded on them.		
Sun's	–	3 to 4 Sun workstations are available for data workup in rms. 2207 and 2219A.		
	–	NARN, VORLON, and XIRTH are hosts for the U500, I500 and I600 instruments, respectively.		
	–	VNMR 6.1c, Felix95, NMRView, NMRPipe, and XEasy software are available on the Sun workstations.		
SGIs	–	NATOTH (Avance host computer) and GQUAN (for off-line data workup)		

II. Facility Personnel

Director, Chemistry Instrument Center

Prof. Paul M. Treichel Rm. 6359A 262-8828 treichel@chem.wisc.edu

Director, Magnetic Resonance Facility

Dr. Charles G. Fry Rm. 2128 262-3182 o fry@chem.wisc.edu
(Charlie) 276-0100 h cgfry@facstaff.wisc.edu

Associate Director, MRF

Monika Ivancic Rm. 2132 262-7536 o

NMR Engineer

Marv Kontney Rm. 2210 262-0563 marv@chem.wisc.edu

ESR Engineer

Roger Clausen Rm. 2225 262-8196 clausen@chem.wisc.edu

Teaching Assistants

Brian Popp Rm 6128 265-8193 bvpoppp@wisc.edu
Mike Birkeland Rm 2201B 262-7536 dalemowrey@hotmail.com